


## How-to-collect data directly on the server (SMART 1000/FRAMBO/GADDS)

### 1. Start SMART / New Project

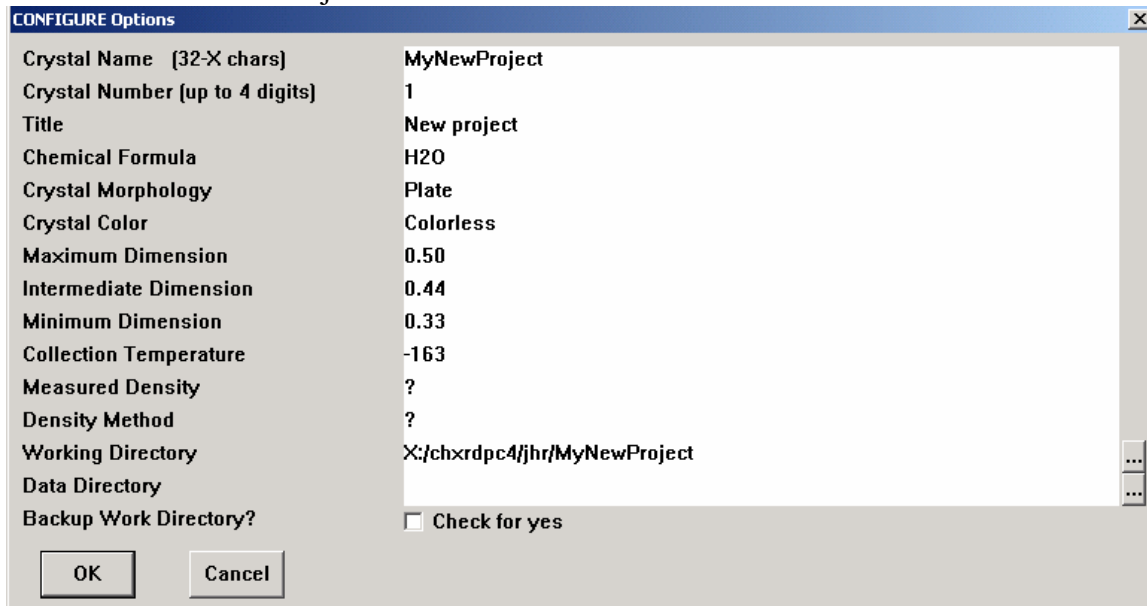

2.

For the working directory type X:/computer\_name/PI\_initials/your\_project\_name

Where computer\_name = chxrdpc4 (SMART1) or chxrdpc5 (SMART2) etc..

PI\_initials = myd, djd, jhr, fpg etc...

your\_project\_name = project

3. Collect data as normal
4. In this case the data is written to  
xray.chem.tamu.edu/data/chxrdpc4/jhr/MyNewProject
5. To reduce the data use SAINT as normal but now input the data from drive X:
6. You may login to xray.chem.tamu.edu and view your data or transfer your data to a new computer.
7. See Lab Managers for User Name and Password on the xray.chem.tamu.edu