

## Use of Radiation Surveillance Equipment.

### Audio "pancake" Geiger Counter.

- 1) Detach detector and remove plastic cover to expose window.
- 2) Turn selection switch to Battery and read status of power. If low replace batteries
- 3) Turn selection switch to lowest detection level.
- 4) Place detector near area to exam and read meter.

### Non-audio Geiger Counter.

- 1) Detach detector and turn detector core to expose window.
- 2) Turn selection switch to Battery and read status of power. If low replace batteries
- 3) Turn selection switch to lowest detection level.
- 4) Place detector near area to exam and read meter.

### Audio Geiger Counter. Small black self-contained device.

- 1) Depress button on front of device. A red light will be seen.
- 2) Place device near area to exam. Read meter and listen to clicks.